

Primary methods for detecting cytosine modifications

Method	DNA prep	Notes
WGBS-Seq Whole Genome Bisulfite-Seq	Na-bisulfite converts non-methylated cytosines to U (U read as T). 5mC and 5hmC are protected from conversion and read as C. Sequencing libraries are prepared by random-primer extension & sequenced with 2x75-100 bp reads lengths.	> 38 million CpG's queried in human. Requires > 90Gb sequence data/sample. Cost ≈ US\$5,000/sample*.
RRBS-Seq Reduced representation Bisulfite seq.	DNA digested with <i>MspI</i> (C CGG), ≈100-150 bp fragments isolated, treated with bisulfite, libraries prepared and sequenced with 1x75 bp.	Queries ≈ 85% CpG islands (≈ 2M CpG's) & 60% of RefSeq promoters. Requires ≈ 40-50M reads, 3-5 Gb/sample. Cost ≈ US\$300/sample*.
MeDIP Methylated DNA immuno-precipitation	DNA sonicated to 100-300 bp, end-repaired, ligated to adapters, and denatured. 5mC containing fragments are captured with anti-5mC magnetic beads, and sequenced (1x75-100 bp). MeDIP may be combined with bisulfite conversion.	5mC detected at ≈150 bp resolution. Bias towards hypermethylated regions. Requires ≈ 60M reads, ≈ 5Gb/sample. Cost ≈ US\$300/sample*.
MIRA Methylated CpG island recovery assay	DNA sonicated to 100-300 bp, end-repaired, ligated to adapters and fragments captured with GST-labeled MBD2b and His-tagged MBD3L1 proteins. Purified by glutathione-magnetic beads. Sequenced with 1x75 bp reads.	Detects mCpGs at ≈ 150 bp resolution. Bias towards hypermethylated CpGs. Non-CpG methylation not detected. Requires ≈ 60M reads, ≈ 5Gb/sample. Cost ≈ US\$300/sample*.
HM450 array Human methylation 450k array	Bisulfite conversion of DNA. C to T changes at defined genomic positions detected by the Infinium assay.	Samples 96% CpG islands, 485,000 CpGs, 99% RefSeq promoters, 3' & 5' UTR, 1st exon, gene body, 3' UTR, shores and shelves. Cost ≈ US\$240/sample*.
Tab-Seq Tet-assisted bisulfite seq	β-glucosyltransferase adds glucose to 5hmC but not 5mC. <i>Tet</i> enzymes oxidize 5mC to 5caC which converts to U on bisulfite treatment but glucosylated 5hmC cannot be oxidized. hmC will be read as C and mC will be read as T.	Detects genome-wide 5hmC at single base resolution. Requires > 90Gb/sample. Cost ≈ US\$5,000/sample*.
hMe-DIP Hydroxy Methylated DNA Immuno Precipitation	DNA sonicated to 100-300 bp, end repaired & ligated to adapters. hmC fragments captured with anti-5hmC labeled magnetic beads. Bound fragments sequenced with 100 bp reads.	Detects hydroxymethylated cytosines at about 150 bp resolution. Requires about 5Gb sequence data/sample. Cost ≈ US\$300/sample*.
RRHP-Seq Reduced Representation Hydroxymethyl-cytosine Profiling Sequencing	DNA digested with <i>MspI</i> (C CGG) & ligated to adapters. hmC glucosylated with β-glucosyltransferase making it resistant to re-digestion with <i>MspI</i> . Treating the library with <i>MspI</i> selectively digests all non-hmC fragments.	≈ 85% of CpG islands (2 M CpG's) & 60% of RefSeq promoter regions queried. Requires ≈ 30-50M reads, 3-5 Gb/sample. Cost ≈ \$300/sample*.

* Cost is provided for guidance only & may not reflect your actual cost.

Techniques in brief

COBRA. Combined Bisulfite followed by Restriction Analysis. RE recognition affected by bisulfite treatment.

C-Subtraction. *MseI* (methylation independent) digestion, ligation to linkers, then digestion with methylation sensitive *BstUI* or *HpaII* to reduce unmethylated DNA.

HELP. HpaII enrichment by ligation PCR. Differential restriction of methylated and unmethylated CpG sites.

MCA. Methylated CpG-island amplification. DNA cut with *SmaI* (blunt ends, cannot cut mC) followed with *XmaI* (sticky ends, cuts mC and C). Sticky end adapters select for mC fragments.

MeDIP. Methylated DNA immunoprecipitation with anti-methylcytosine antibody.

Methyl light. Methylation specific qPCR (see MSP).

Methylation trapping of methyl transferases onto DNA with 5-Aza-2'-deoxycytidine (decitabine) followed by immunoprecipitation.

MIRA. Methylated CpG island recovery assay. Capture with MBD2b and MBD3L1 protein heterodimer.

MSDK. Methylation specific digital karyotyping. Cleavage with methylation sensitive *AscI*. Sequence tags sequenced and mapped like SAGE.

MSP. Methylation specific PCR on bisulfite converted DNA with two sets of primers to amplify either -C or -T base.

OxBS. K-perruthenate oxidizes only 5hmC to 5fC that is converted to U by bisulfite.

RLGS. Restriction Landmark Genome Scanning with 2D gel of methylation sensitive *NotI* and *AscI* restricted DNA.

RRHP. Reduced Representation Hydroxymethylation Profiling.

RRBS. Reduced Representation Bisulfite Sequencing.

Tab-Seq. Tet-assisted bisulfite sequencing of 5-hmC.

WGBS-Seq. Na-bisulfite conversion of C → T. mC and hmC are not converted and are read as C by whole genome next generation sequencing.

References:

- Harris RA et al (2010) Comparison of sequencing-based methods to profile DNA methylation and identification of monoallelic epigenetic modifications. *Nature Biotech* 28, 1097-1105.
- Guibert S, Weber M (2013) Functions of DNA methylation and hydroxymethylation in mammalian development. *Curr Top Dev Biol* 104, 47-83.
- Laird PW (2010) Principles and challenges of genome-wide DNA methylation analysis. *Nature Rev Genetics* 11, 191-203.
- Meissner A (2012) Epigenomics, special issue. *Genome Biol* 13(10).
- Stevens M et al (2013) Estimating absolute methylation levels at single CpG resolution from methylation enrichment and restriction enzyme sequencing methods. *Genome Res.* 10.1101/gr.152231.112.
- Xie W et al (2013) Epigenomic analysis of multilineage differentiation of human embryonic stem cells. *Cell* 153, 1134-48.
- Yu P (2013) Spatiotemporal clustering of the epigenome reveals rules of dynamic gene regulation. *Genome Res* 23, 352-64.
- Yu M et al (2012) Tet-assisted bisulfite sequencing of 5-hydroxymethylcytosine. *Nature Protocols* 7, 2159-70.

FOR RESEARCH USE ONLY

© 2013 Illumina, Inc. All rights reserved.

Illumina, IlluminaDx, BaseSpace, BeadArray, BeadXpress, cBot, CSPPro, DASL, DesignStudio, Eco, GAllx, Genetic Energy, Genome Analyzer, GenomeStudio, GoldenGate, HiScan, HiSeq, Infinium, iSelect, MiSeq, Nextera, NuPCR, SeqMonitor, Solexa, TruSeq, TruSight, VeraCode, the pumpkin orange color, and the Genetic Energy streaming bases design are trademarks or registered trademarks of Illumina, Inc. All other brands and names contained herein are the property of their respective owners.